


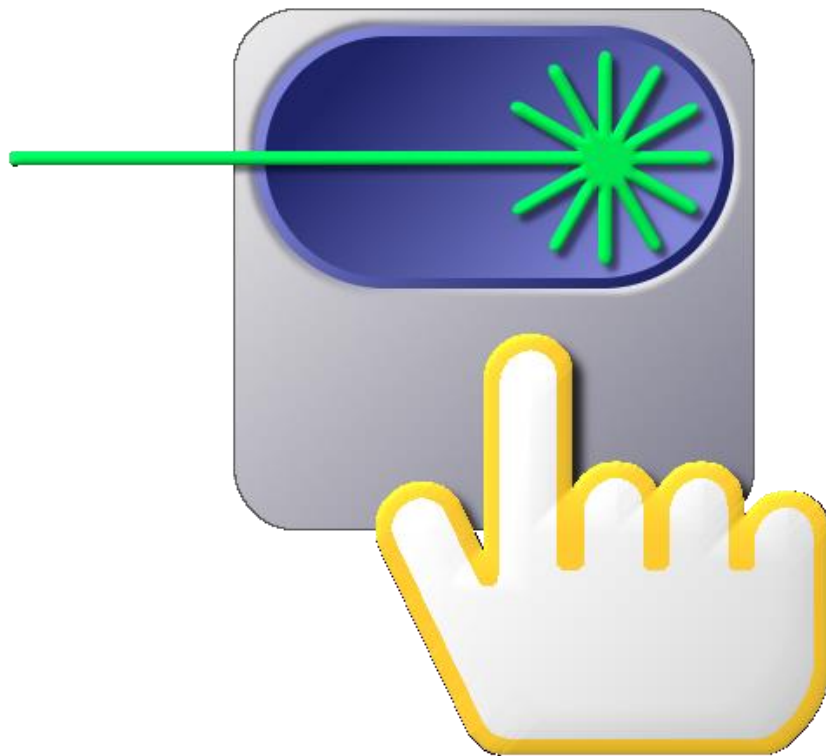

# **Laser Control Software (LabView)**

## **User Manual**



**EdgeWave GmbH**  
Innovative Laser Solutions

# EdgeWave Laser Control Software (LabView)

Copyright c 2010 EdgeWave GmbH  
Doc. Ver. 1.0.3 - February 18, 2017

Every effort has been made to ensure that the data given in this document is accurate. The information, figures, tables, specifications and schematics contained here in are subject to change without notice. EdgeWave makes no warranty or representation, either expressed or implied with respect to this document. In no event will EdgeWave be liable for any direct, indirect, special, incidental or consequential damages resulting from any defects in its documentation.

This manual is based on EdgeWave Laser Control version 1.1.0 and above. The version of the software can be found in the settings tab by clicking on the info button.

EdgeWave GmbH  
Carlo-Schmid-Str. 19  
D-52146 Würselen  
Germany  
Tel: +49(0)2405-4186 0  
Fax: +49(0)2405-4186 33  
Email: [info@edge-wave.com](mailto:info@edge-wave.com)  
Internet: <http://edge-wave.com>

# Table of Content

1. Safety.....	1
2. Installation.....	2
2.1 File list .....	2
2.2 Requirement.....	2
2.2.1 Hardware requirement .....	2
2.2.2 Software requirement.....	2
2.3 Software installation and uninstalling .....	3
2.3.1 Installation .....	3
2.3.2 Uninstalling .....	3
3 Operation .....	4
3.1 Main window .....	4
3.2 Terms .....	8
3.3 Regular Operation .....	14
3.3.1 Laser Operation Panel .....	14
3.3.1.1 Laser Operation Panel (Nano Second Laser).....	14
3.3.1.2 Laser Operation Panel (Pico Second Laser) .....	17
3.3.2 Tab Control Panel .....	20
3.3.2.1 Current Control Tab.....	20
3.3.2.2 Oven Control Tab .....	21
3.3.2.3 Advanced Info Tab .....	22
3.3.2.4 Settings Tab.....	24
3.3.2.5 Chiller Control Tab.....	25
4 Options.....	28
4.1 Laser Connection Settings.....	28
4.2 Hot Keys .....	30
4.3 Preset Currents .....	30
4.4 Save Settings.....	31



### **1. Safety**

The EdgeWave Laser Control Software is intended to control an EdgeWave INNOSLAB laser through a serial (RS232) port of the computer. Laser radiation may affect a person's health or may otherwise cause damage. Prior to installation and operation compliance with all relevant laser safety regulation has to be secured. The client shall solely be responsible to strictly comply with all applicable and relevant safety regulations regarding installation and operation of the system at any time.

## 2. Installation

### 2.1 File list

The EdgeWave Laser Control Software is delivered with the following file(s):

- Laser Control.exe

Laser Control which provides the full control to the Laser via RS232 (serial) port.

- Settings.ini

The file used to save all settings for the EdgeWave Laser Control Software.

- Data (Folder)

LabView runtime engine needs the DLL-files in this folder to run the Laser Control Software.

### 2.2 Requirement

#### 2.2.1 Hardware requirement

- The EdgeWave Laser Control Software is designed to control EdgeWave INNOSLAB Laser. An EdgeWave Laser with essential peripheral equipment, e.g. EdgeWave Power Supply, chiller, trigger generator, etc. is required.



Please read the manual of EdgeWave Laser carefully and connect all cables properly prior to use EdgeWave Laser Control Software.

- A computer to control the Laser.
  - A serial (RS232) port is required to control an EdgeWave Laser with EdgeWave Laser Control Software. Both on-board com port and USB to RS232 adapter works.
  - 10MB free hard drive space.
  - The software itself will use less than 50MB memory when running, which can be provided by any modern computer.
  - A monitor with a resolution of min. 1280x720 pixels.

#### 2.2.2 Software requirement

- EdgeWave Laser Control Software works only in Windows™ platform, including both 32 and 64 bits version of Windows XP, Windows Vista, Windows 7 or above.
- LabView run-time engine 2010 (32-bit standard RTE) and NI-Visa 5.1 RTE or above. (The RTE can be downloaded from Ni homepage [www.ni.com](http://www.ni.com))

- Windows user privilege
  - The EdgeWave Laser Control Software writes all settings into *Settings.ini* in the same directory when it is closed. The setting file will be created automatically in case of missing. The Modify, Create and Read privilege to file *Settings.ini* must be granted to current user to avoid some warning message.
  - The software controls an EdgeWave Laser via the serial (RS232) port. The read and write privilege to the using serial port has to be granted.

### 2.3 Software installation and uninstalling

The EdgeWave Laser Control Software is a green software which means NOTHING will be written into the registry or system folder by itself.

#### 2.3.1 Installation

1. Copy all files delivered into the destination directory.
2. Create a shortcut to the preferred location. For instance, a shortcut will be created on the desktop of Windows by right clicking on *Laser Control.exe* with the mouse and then select "Send To" → "Desktop (create shortcut)", as shown below:

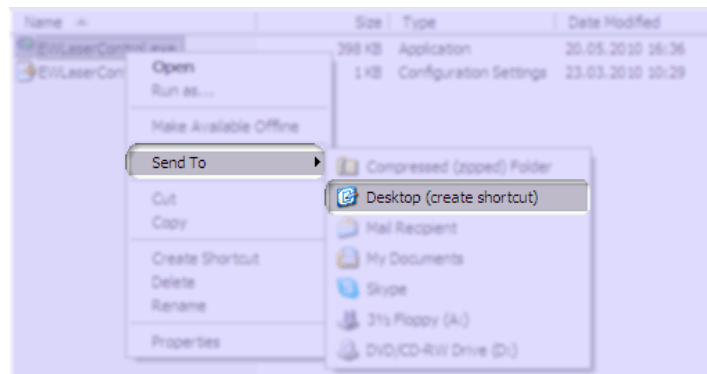


Figure 1: Create a shortcut on desktop

#### 2.3.2 Uninstalling

1. Delete the shortcut pointed to *Laser Control.exe*.
2. Delete *Laser Control.exe*, *Settings.ini* and any other files and folders copied with this software under the directory of EdgeWave Laser Control.

## 3 Operation

### 3.1 Main window

When launch the Laser Control Software for the first time, it will try to connect EdgeWave Laser via a serial port. If no EdgeWave Laser is connected (Figure 2), it will try to reconnect automatically for every 5 seconds If *search for laser periodically* option is activated (checked). Both serial port and the auto-connect feature are adjustable in the *Settings* tab.

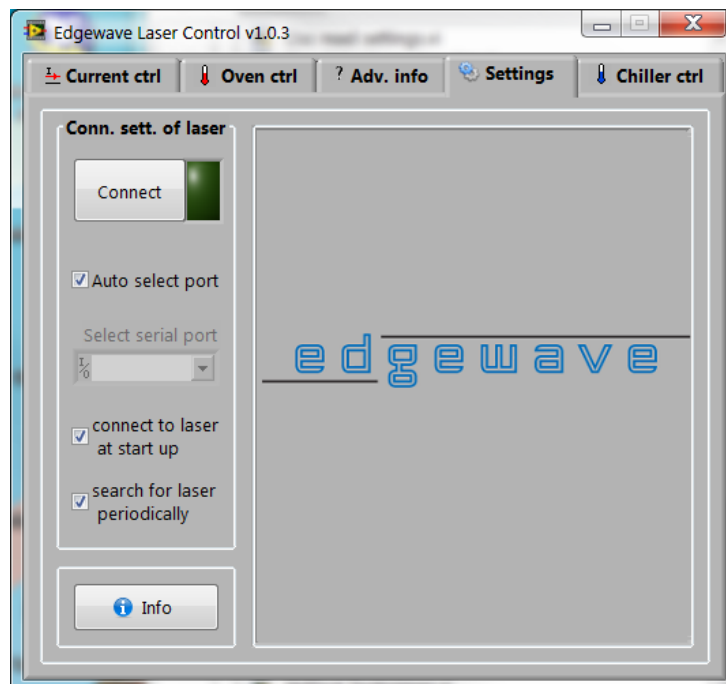


Figure 2: Settings tab – Laser disconnected

To establish a connection to the laser, the following requirement must be fulfilled:

- An EdgeWave laser is connected to a working serial port of the computer.
- The RS232 cable used to connect the laser and the computer must be a 1-to-1 cable.
- Serial port number has to be specified in the laser control software correctly. If the *Auto select port* option is not checked.
- Emergency stop button of the laser has to be released.
- Safety loops on the rear panel of laser power supply is connected properly and the corresponding switches are closed. Please find more details about safety loops in the user manual of EdgeWave Laser.
- The Laser power supply is connected to the mains and switched on.



### 3. Operation

Once the laser is found and connected, *Laser Operation* panel (Figure 3) will appear. This panel provide an easy way to set most of laser parameters and displays warnings or errors. The Current Control tab (Figure 3) displays the current status. Please refer to section 3.3.1 for detailed information.

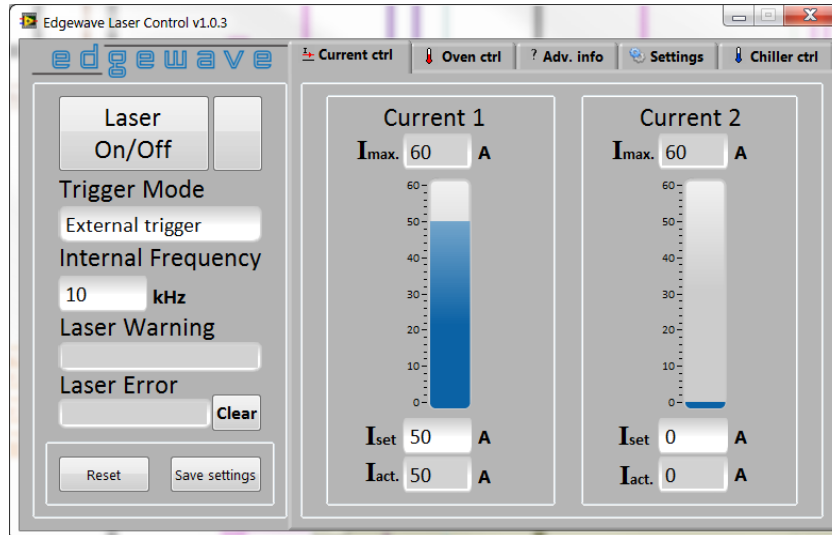


Figure 3: Laser operation panel and current control tab – Nanosecond laser

If a pico second laser is connected, advanced setting for pico second laser (Figure 4) will appear. The high voltage level of the modulator, modulator mode and the pulse picker settings can not be changed while the laser is on. Please refer section 3.3.1.2 for more information.

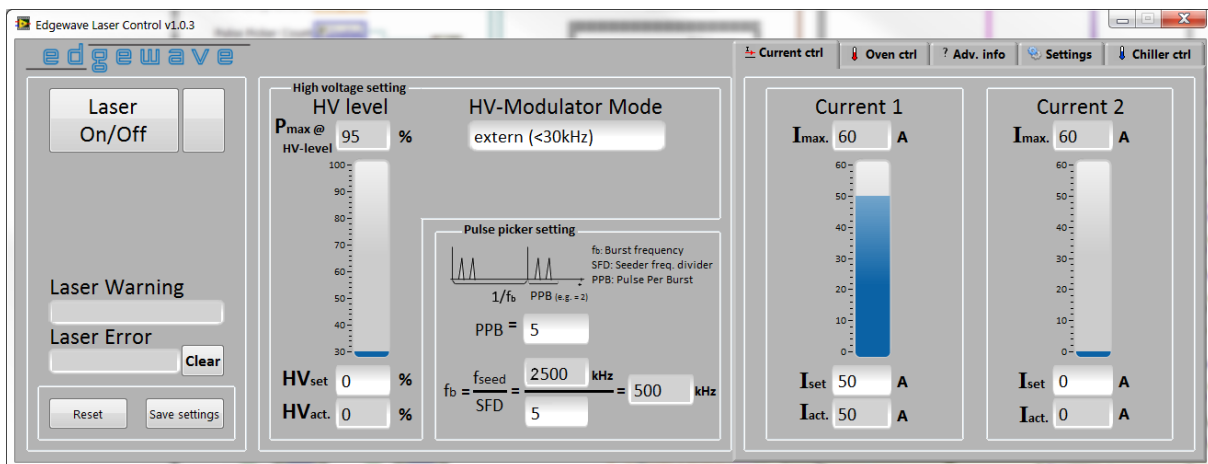


Figure 4: Laser operation panel and current control tab – Pico second laser

### 3. Operation

In the Oven Control tab (Figure 5) the activated ovens will be shown. Ovens will be activated automatically during connecting the laser. In the Advanced Info tab (Figure 6) there is the possibility to activate or deactivate the ovens individually. Please refer to section 3.3.2.2 for more information.

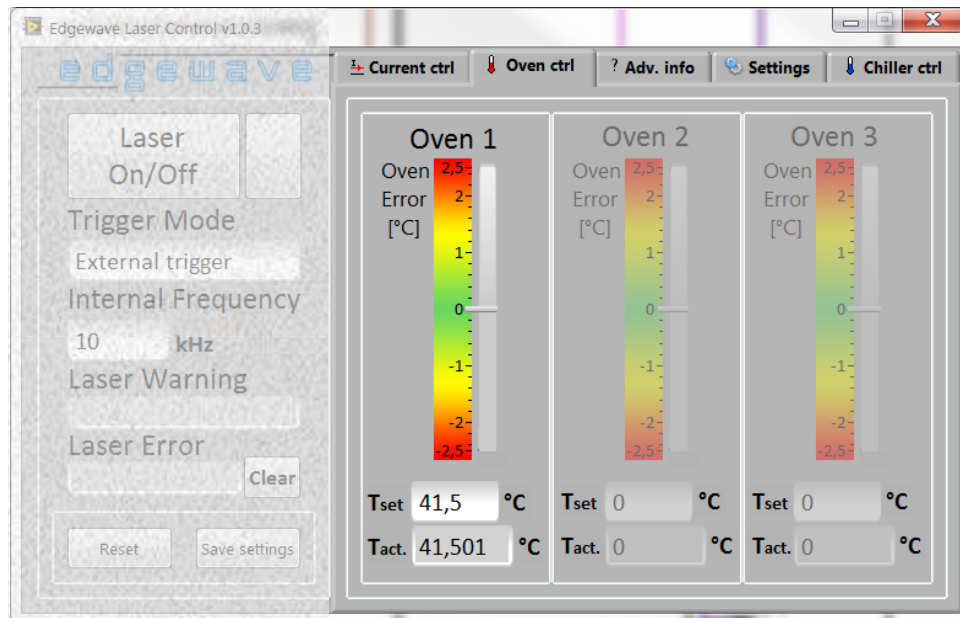


Figure 5: Oven control tab

Information (e. g. Serial number, humidity or cooling water settings) about the connected laser are displayed in the Advanced Info tab (Figure 5). In this tab a terminal can be activated, with which all control commands of EdgeWave Laser can be executed. Please refer section 3.3.2.3.

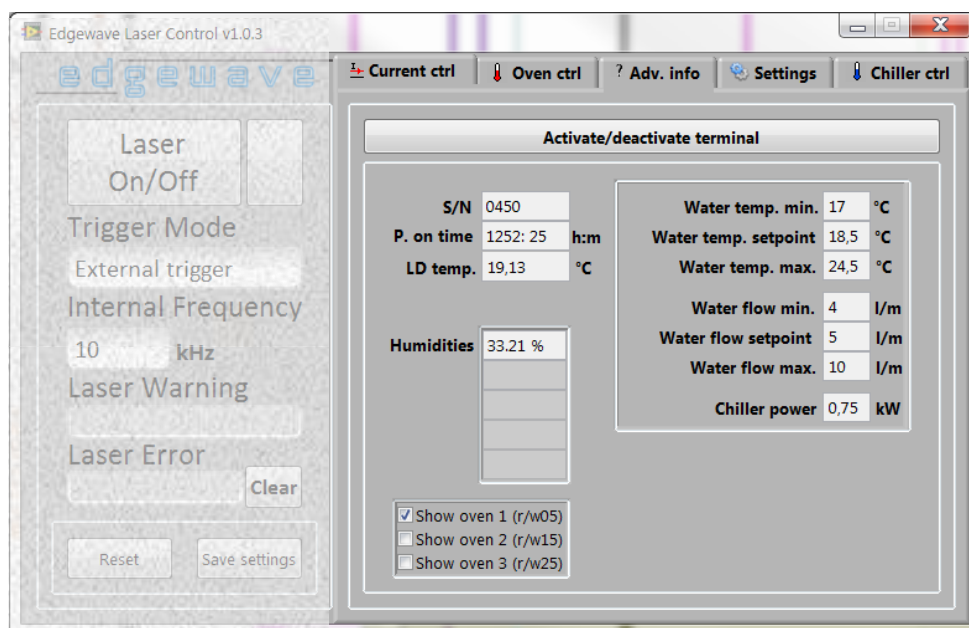


Figure 6: Advanced info tab – Nanosecond laser

A Termotek chiller with a serial interface (RS232) can be controlled over the *Chiller Control* tab (Figure 7). It allows a simple control of a chiller (e. g. water temperature set point and turn on/off the chiller) and displays the outlet water temperature and water flow rate.

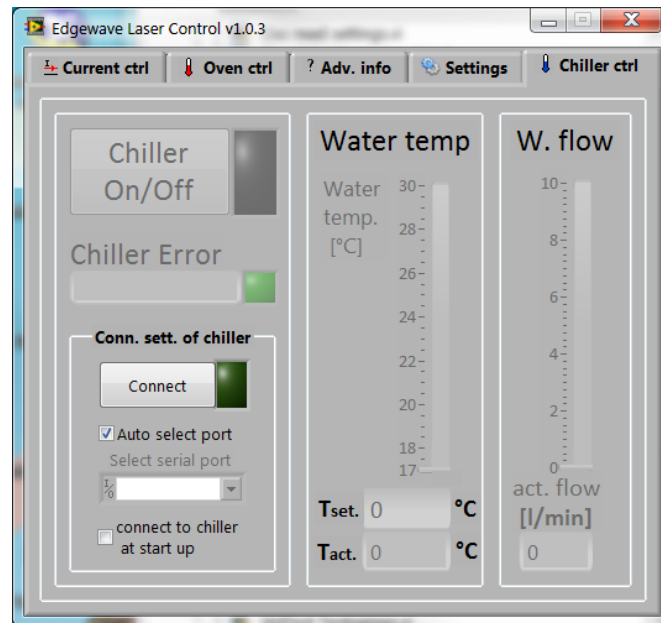


Figure 7: Chiller control tab. – chiller disconnected

To establish a connection to a chiller, the following requirement must be fulfilled:

- A Termotek chiller is connected to a working serial port of the computer.
- The RS232 cable used to connect the laser and the computer must be a 1-to-1 cable.
- Serial port number has to be specified in the chiller control tab correctly. If the *Auto select port* option is not checked.
- The chiller is connected to the mains and switched on.

### 3.2 Terms

The terms below are subject to the use of EdgeWave Laser. Their definition is important in understanding the operation procedure, trouble shooting and so on. Some of the definitions are valid only inside this manual and during operating EdgeWave Laser.

#### Laser output

Laser output means required laser beam emitting from the output aperture of the laser head. To have laser output, the following must be fulfilled if the other necessary parts (e.g. cooling water) are configured properly and working well:

1. Laser ON
2. The current is higher than the threshold
3. Proper TTL trigger signal is provided if working in External trigger mode
4. Proper TTL trigger signal for external modulator is provided if required

#### Laser ON/Off

When Laser On/Off is set, the control board has to prepare some initializing procedures for laser output, e.g. set the current of Laser diodes to nominal value and turn the shutter to output position if present.



The diode current and pulse picker settings of a pico second laser can not be changed and will be deactivated while laser on.



Laser ON should be considered more like "Laser is ready for ON". To have laser beam output, all conditions for Laser output described above must be fulfilled.

When laser is Off, current of Laser diodes is set to 0A despite of the nominal current and the shutter is turned to off position (if present).

#### Q-Switching

Most of EdgeWave lasers are working at Q-switching mode, which is a technique to produce pulsed output beam with high peak power and short pulse width. It is achieved by putting an optical switch inside the laser's optical resonator. Normally there is not pulse output when the switch is off, while a giant pulse is produced once it switches to on.

#### Trigger

The TTL signal used to control the on/off state of the Q-switch. It controls the repetition rate and the timing of laser pulses. It also regulates the pulse energy in some kinds of EdgeWave laser.

### Trigger mode

Control mode of the Q-switch in EdgeWave laser. There are 3 kinds of mode: CW (continuous wave), External trigger and Internal trigger.

#### External trigger

The trigger signal is generated outside of EdgeWave laser and coupled into the laser through the trigger port of the laser. This is the most favourable mode for most applications because the user can control the timing and the other features.

#### Internal trigger

The trigger signal is generated by the control board of EdgeWave laser. It is not controllable by user. For some model the internal trigger signal might becomes unstable at >5kHz.

#### CW

The Q-switch does not work under CW mode. The laser resonator will have constant output if there is enough pump power.

### Internal Frequency

The internal trigger frequency is generated by the control board in laser.



In some lasers equipped with certain first pulse killer, the frequency set here also tells the controller to perform the pulse optimizing according to this frequency. For such laser, the maximum working frequency must be set here. Otherwise the laser might be locked down at lower power level, or pulse energy is unstable.

### Laser status indicator

The status indicator on the right side of the *Laser On/Off* control displays the actual status of the laser analogue to the status LED on the laser head and power supply of some lasers.



The blue color of the *Status indicator* displays that the laser is ready to use.



If the laser is on the color of the *Status indicator* changes to green



On any error the *Status indicator* changes its color to red.

### High voltage settings (Pico second laser)

The high voltage level of the modulator regulates the avg. laser power of pico second laser. While the laser is on, the set point of the current can not be changed. By changing the high voltage level from 30% to 100%, the avg. power of the laser can be manipulated.



Please note that the high voltage level does not correspond with the avg. laser power. The *Pmax@HV-level* displays the high voltage level set point at which max. laser power can be emitted.

### HV-Modulator mode (Pico second laser)

Control mode of the modulator in EdgeWave pico second laser. There are 4 kinds of mode: HV Always on, HV Always off, External and Internal.

#### Internal

The internal signal is generated by the control board of laser. The maximum frequency is displayed on the *HV-Modulator Mode* control element. If changed to Internal mode an input field for the duty cycle frequency will appear.

#### External

The signal is generated outside of EdgeWave laser and coupled into the laser through the trigger port of the laser. This is the most favourable mode for most applications because the user can control the timing and the other features. An external function generator can be synchronized through the sync. output of the laser head. The maximum frequency is displayed on the *HV-Modulator Mode* control element.

#### HV Always off

The modulator is turned off. No pulses will be emitted. This is the best way to turn off the laser power for a short time.

#### HV Always on

The modulator is on. This means that every laser pulse will be emitted which is picked by the pulse picker.

#### Duty cycle frequency

The duty cycle signal can be generated by the laser control board or by an external function generator. Laser pulses will emit only in the positive width of the signal. The maximum frequency, which can be used, is displayed on the *HV-Modulator Mode* control element.



The duty cycle signal can be used only if HV-modulator mode is switched to internal or external mode. If modulator mode changed to internal mode an input field for the duty cycle frequency will appear.

#### Pulse picker settings (Pico second laser)

For different applications it is possible to set a repetition rate lower than the seeder frequency. An AOM (acousto-optic modulator) allows to pick out one or more pulses from a train of pulses (Figure 8) generated by the seeder.

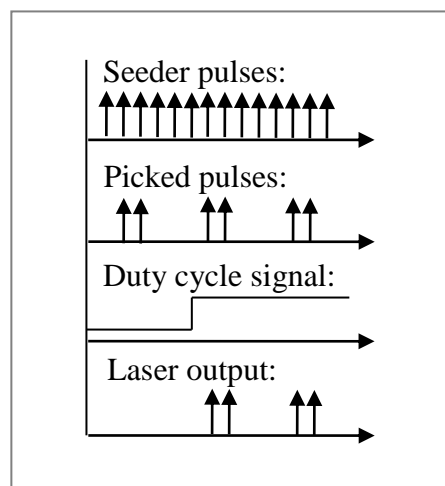


Figure 8: Pulse picker and modulator states

#### Seeder frequency ( $f_s$ )

The *seeder frequency* is displayed on the *pulse picker settings* panel. It helps to calculate the *burst frequency*.

#### Burst frequency ( $f_b$ )

The *burst frequency* is calculated by dividing the *seeder frequency* with the *seeder frequency divider*. Every burst can consist of one or more pulses. The pulse count can be set over *pulse per burst (PPB)* input field.

**Seeder frequency divider (SFD)**

An integer value can be set to divide the seeder frequency. To set a different SFD laser must be off.



Please note that by higher SFD values the amplifier current (I2) is limited! To be able to set a higher SFD the amplifier current must be set correct. A frequency current table on the Advanced Info tab (Figure 9) shows the possible settings.

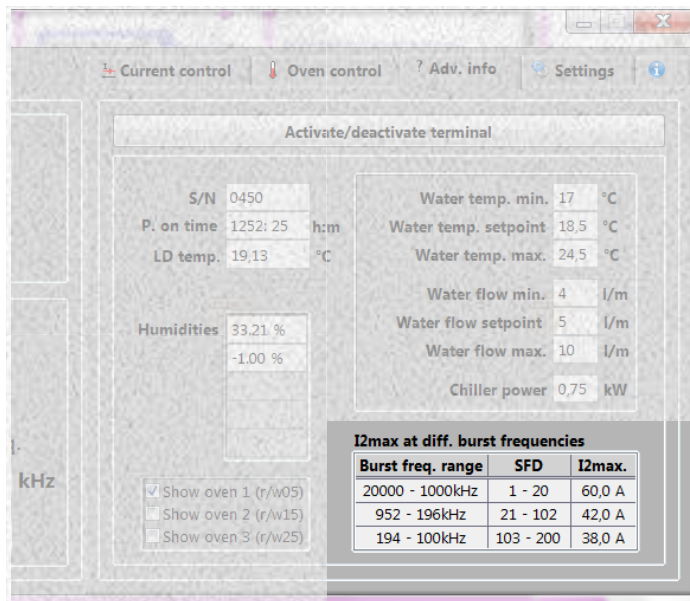


Figure 9: Frequency current table – Advanced Info tab

**Pulse per burst (PPB)**

An integer value can be set to change the pulse count per burst. If PPB is zero no pulse will emit!



#### Hyper Terminal

A windows application for terminal emulation which supports communication over TCP/IP networks, Dial-Up Modems, and serial ports (RS232). It was bundled with Windows from version Windows 95 to Windows XP.

This tool is useful for troubleshooting. There are also some other software provides similar functions.

The function of hyper terminal to communicate with a connected EdgeWave laser can be used by activating the terminal in the Advanced Info tab (Figure 10). Please refer section 3.3.2.3 for information about using the terminal function.

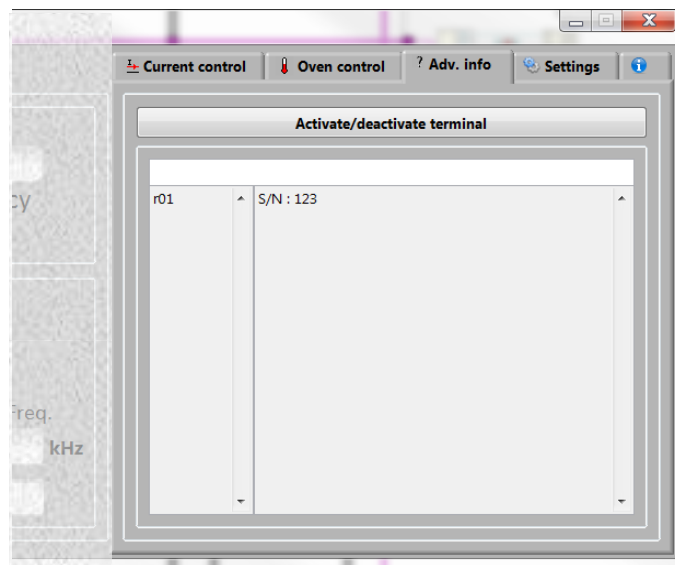


Figure 10: Advanced info tab with activated terminal

#### Info

Displays the version number (Figure 11) of the EdgeWave laser control software and contact information.

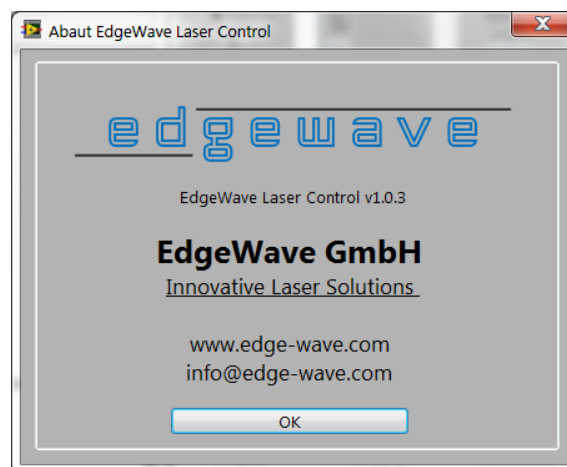


Figure 11: Info dialog

### 3.3 Regular Operation

The regular operation window provides a GUI (graphical user interface) to interact with an EdgeWave laser. Only a few value inputting and mouse clicking are required with such panel to control the Laser. Most of laser operation can be done here.

#### 3.3.1 Laser Operation Panel

*Laser Operation* panel is located at the left side of the window. If a nano second laser is connected (Figure 12), the panel will be visible. If a pico second laser is connected, advanced controls and indicators will appear (Section 3.3.1.2).

##### 3.3.1.1 Laser Operation Panel (Nano Second Laser)

The Laser Operation panel for nano second laser consists of the input fields for *Trigger Mode*, *Internal Frequency*, indicators for *Laser Status*, *laser warning* and *error*.

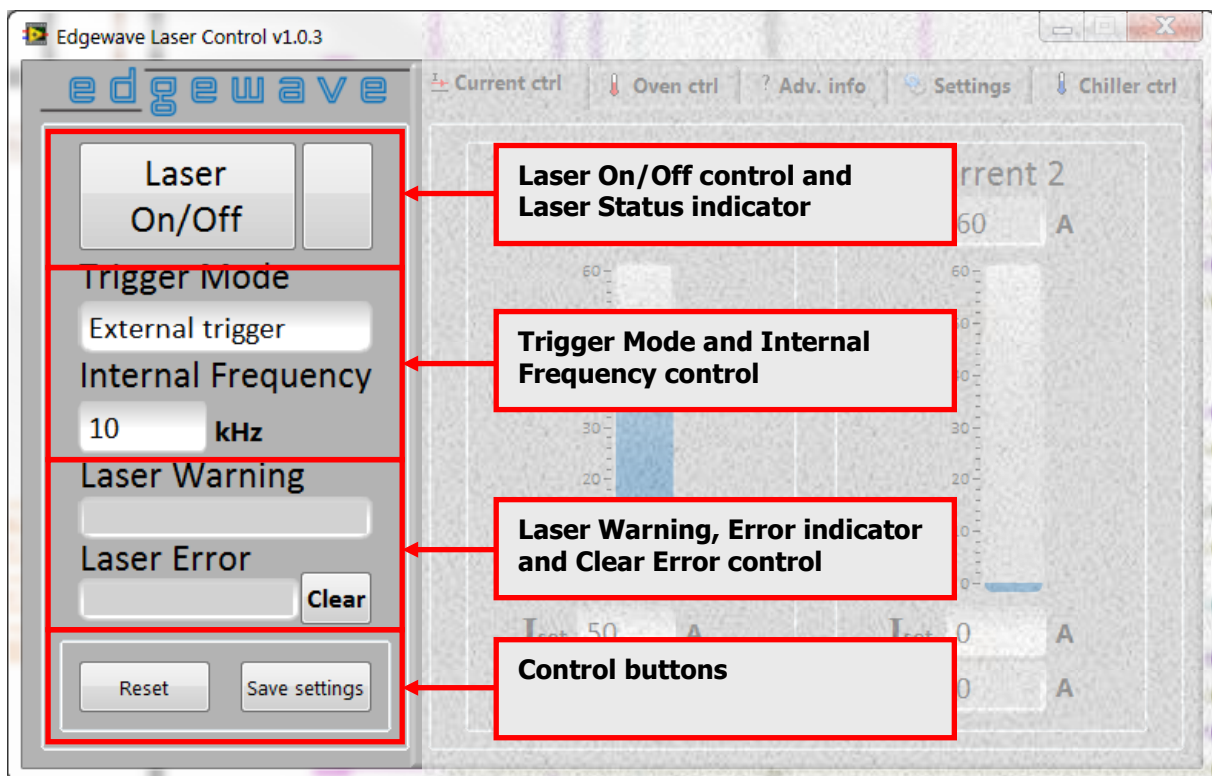


Figure 12: Main operation panel


### Laser On/Off Control

To turn on or off the laser. Please refer to section 3.2 for the definition of laser on and off

### Laser Status Indicator

It shows the laser on, off and error status of the laser. Please refer to section 3.2 for detailed information.

### Trigger Mode Control

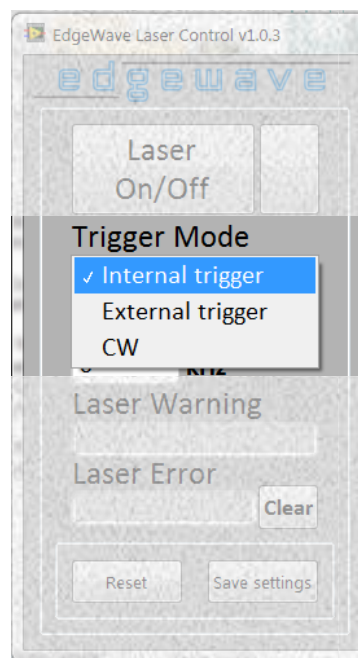

The trigger mode can be changed easily by clicking the corresponding control (Figure 13). The trigger modes are: Internal, external and CW (continuous wave). Please refer to section 3.2 for the definition of trigger modes.

Figure 13: Trigger Mode control

### Internal Frequency

An internal trigger frequency in kHz can be input here.



In some Lasers equipped with certain first pulse killer, the frequency set here also tells the controller to perform the pulse optimizing according to this frequency. For such Laser, the maximum working frequency must be set here. Otherwise the Laser might be locked down at lower power level, or pulse energy is unstable.

### **Laser warning**

It shows laser warning messages. If the warning limits reached (e. g. water temperature too high or low) a warning text will be displayed.

### **Laser Error**

It shows laser error messages. If an error occurs the laser will turn off and *Laser Status* indicator becomes red. Laser is locked down and can not be turn on until the error not solved. To release the laser from lock down state the *Clear error* button must be used.

### **Clear Error**

If laser error happened and solved, click this button to try to clear error message and release the laser from lock down state.

### **Reset**

Click the reset button to perform a cold start. By doing this the laser will be turned off (if it is on) and all settings will be restored to the default value saved before (using *Save Settings* button)

### **Save Settings**

Save all settings (including trigger frequency and mode, nominal current, etc.) into laser controller as default values. Once saved, those settings will be restored after next reset or starting up the power supply.

### 3.3.1.2 Laser Operation Panel (Pico Second Laser)

In case of connecting a pico second laser, advanced controls and indicators will appear (Figure 14).

The pico second laser panel consists of two sub panels:

High voltage modulator panel and pulse picker settings panel.

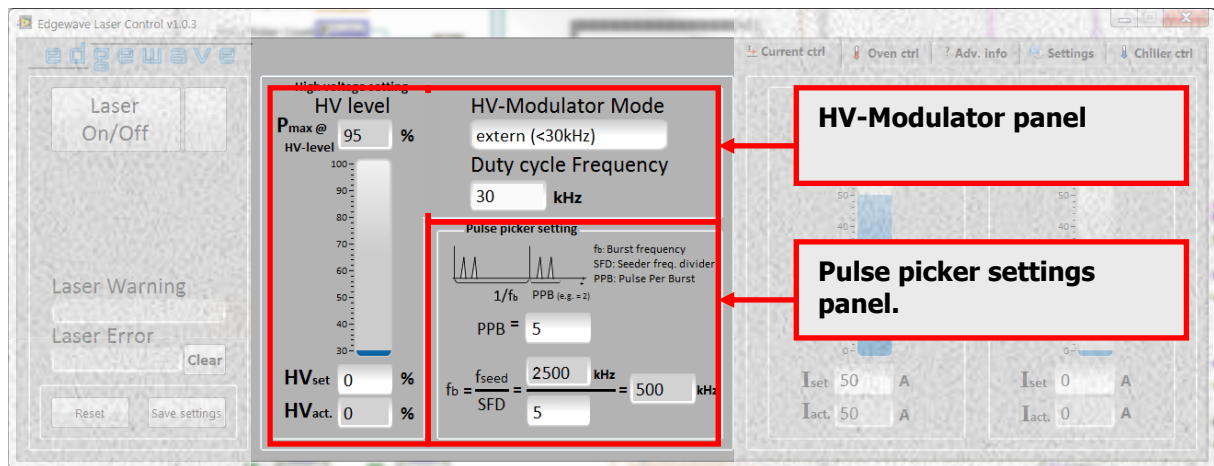


Figure 14: Main operation panel – Pico second laser

#### High Voltage Level Panel

The high voltage level of the modulator regulates the avg. output power of pico second laser. Please refer section 3.2 for more information.

#### Pmax @ HV-level

It shows the high voltage level at which the avg. output power of the laser has its maximum.

#### HVset

High voltage level should be input in this input field.

#### HVact.

It shows the actual high voltage level

## HV-Modulator Mode Panel

### HV-Modulator Mode

Control mode of the modulator in EdgeWave pico second laser. There are 4 kinds of mode (Figure 15): HV Always on, HV Always off, External and Internal. Please refer section 3.2 for detailed information about the modulator modes.

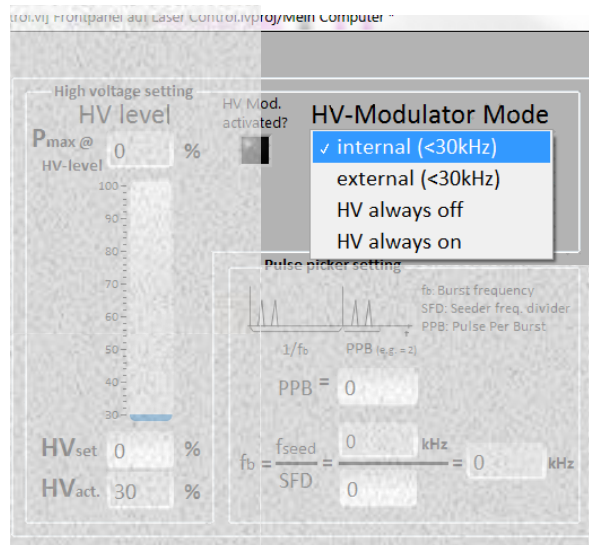


Figure 15: High Voltage Modulator panel

### Duty Cycle Frequency

The duty cycle frequency for the internal modulator mode can be input here. Please refer section 3.2 for more information.

### HV Mod. Activated?

It shows if the high voltage modulator is activated. The modulator can be activated or deactivated with the command:

w178 X (X = 0 to deactivate, X = 1 to activate modulator)



If HV. modulator is deactivated, the laser can not emit any pulses.

### Pulse Picker Settings Panel

This panel allows to set seeder frequency divider and shows the burst frequency (Figure 16). Please refer section 3.2 for the definitions and detailed information.

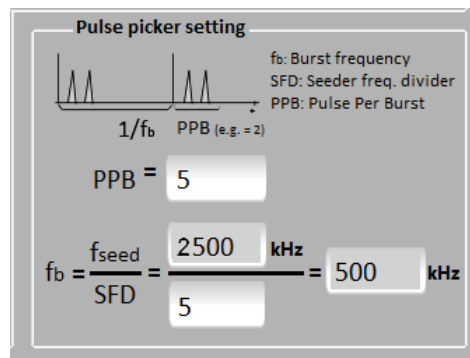


Figure 16: Pulse Picker Settings panel

#### f<sub>seed</sub> (Seeder Frequency)

The seeder frequency is displayed here. It is needed to calculate the burst frequency.

#### f<sub>b</sub> (Burst Frequency)

The burst frequency is calculated by dividing the seeder frequency with the seeder frequency divide (SFD).

#### SFD (Seeder Frequency Divider)

An integer value from 1 to 16 000 can input here. It divides the seeder frequency to allow repetition rates lower than the seeder frequency. Nominal value is sent to the controller only after press enter key in the edit box.



Please note that by higher SFD values the amplifier current (I<sub>2</sub>) is limited! To be able to set a higher SFD the amplifier current must be set correct. A frequency current table on the Advanced Info tab (Figure 9 and 19) shows the possible settings.

#### PPB (Pulse Per Burst)

Every burst can consist of one or more pulses. The pulse count can be set from 0 to SFD value here.



The PPB value must be lower or equal to SFD. If PPB is zero, no pulses will emit!

### 3.3.2 Tab Control Panel

The *tab control panel* consists of five tabs:  
Current Control, Oven Control, Advanced Info, Settings and Chiller Control.

#### 3.3.2.1 Current Control Tab

The current applied on the laser diodes regulates the output power of the optical resonator in a diode pumped laser. For most applications with laser, the power level or pulse energy achieved from laser head has to be controlled through the diode current. The appearance of the *Current Control* tab varies with the hardware configuration of the connected laser (Figure 17).

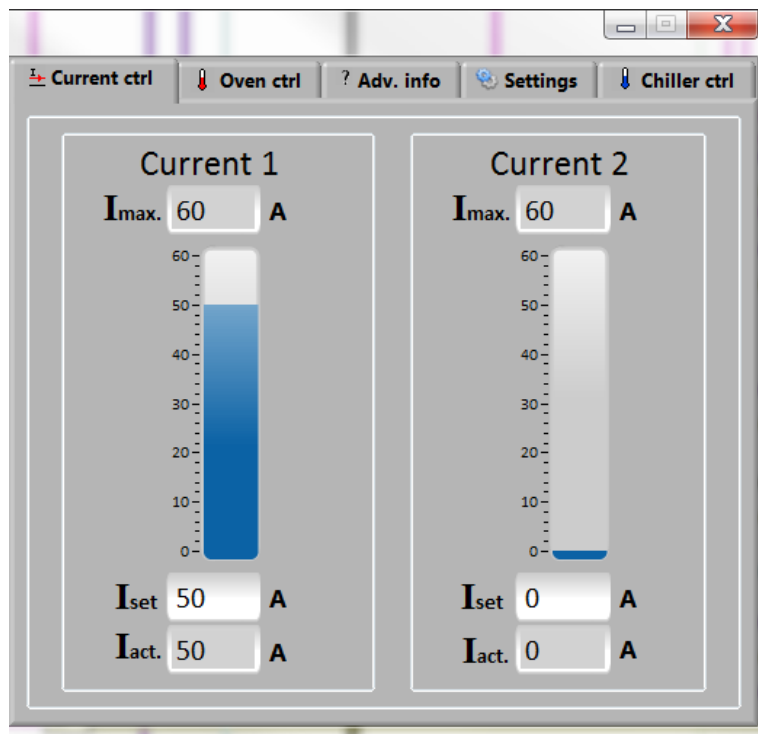


Figure 17: Current control tab – two branches

#### **I<sub>max</sub>**

I<sub>max</sub> (maximum current) shows the maximum possible current settings.

#### **I<sub>set</sub>**

Currents in ampere can be set from 0 to *I<sub>max</sub>* here. Nominal value is sent to the controller only after press Enter key in the edit box.

#### **I<sub>act.</sub>**

A fill slide and an indicator show the actual current in ampere.



### 3.3.2.2 Oven Control Tab

Some EdgeWave lasers have LBO crystal to generate second harmonic of the infrared light (SHG, also known as frequency doubling, e. g. 1064nm -> 532nm). It is also possible to generate laser light with different wavelengths by using more ovens.

The second harmonic generation is very temperature sensitive. To have the best conversion efficiency, it is important to set the optimum oven temperature in which is the LBO crystal embedded.

In the *Oven Control* tab (Figure 18) the activated ovens will shown. Ovens will be activated automatically during connecting the laser. In the Advanced Info tab there is the possibility to activate or deactivate the ovens individually.

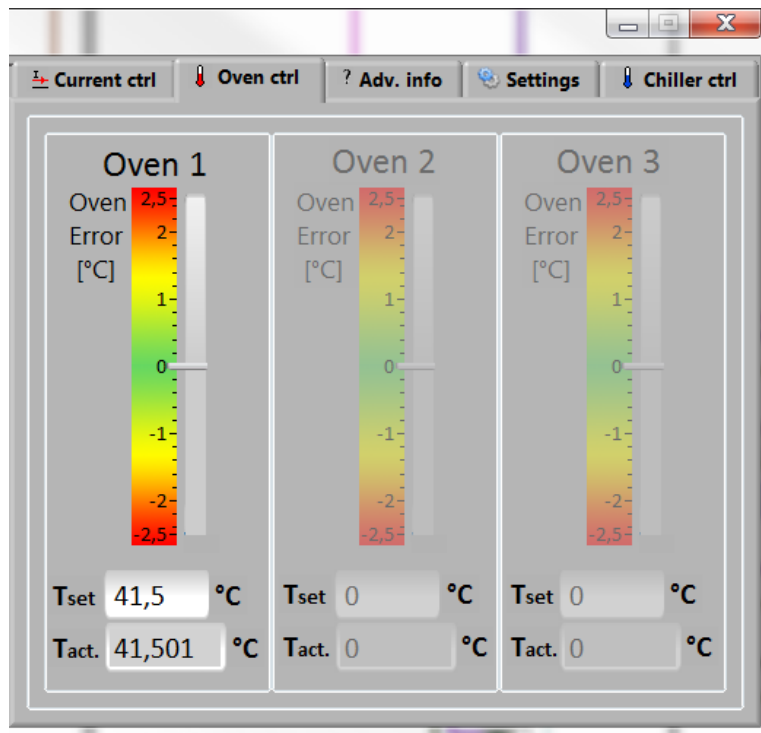


Figure 18: Oven control tab – oven 1 activated

#### Tset

The temperature set point will read from laser while connection and can be changed here. Nominal value is sent to the controller only after press Enter key in the edit box.

#### Tact.

Actual temperature of oven will be shown here.

#### Oven Error

A slider shows the oven temperature error in a range of  $\pm 2.5^{\circ}\text{C}$ .

### 3.3.2.3 Advanced Info Tab

Advanced information about the connected laser displayed in this tab.

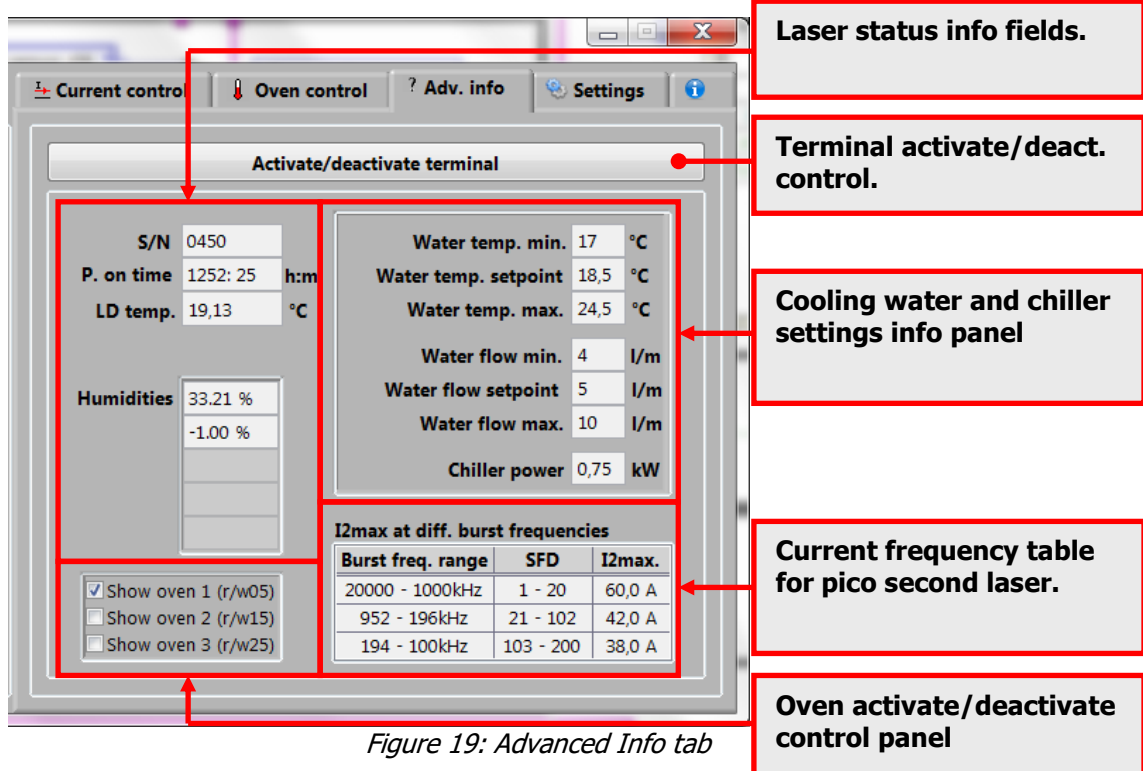


Figure 19: Advanced Info tab

#### Laser Status Info Fields

Laser serial number and different status information will be displayed in these fields.

#### Oven Activate/Deactivate Control Panel

Depended on the hardware configuration of some laser, the auto detection of embedded ovens will not work. In this case one or more ovens can be activated manually here.

#### Cooling Water And Chiller Settings Info Panel

Information about cooling water settings will be shown here. This panel is only visible if it is supported by the connected laser!

#### Current Frequency Table

This panel shows the maximum amplifier current by different SFD settings. Only visible if a pico second laser is connected!

### Terminal Panel

The terminal panel (Figure 20) will be shown by activating it over the *Terminal Activate/Deactivate* control or using hot key (Figure 24). This panel is a terminal-like window for all commands that EdgeWave laser supports. Since the GUI provides basic controls only, the unimplemented commands must be carried out here or use terminal emulation software (e.g. HyperTerminal) instead.

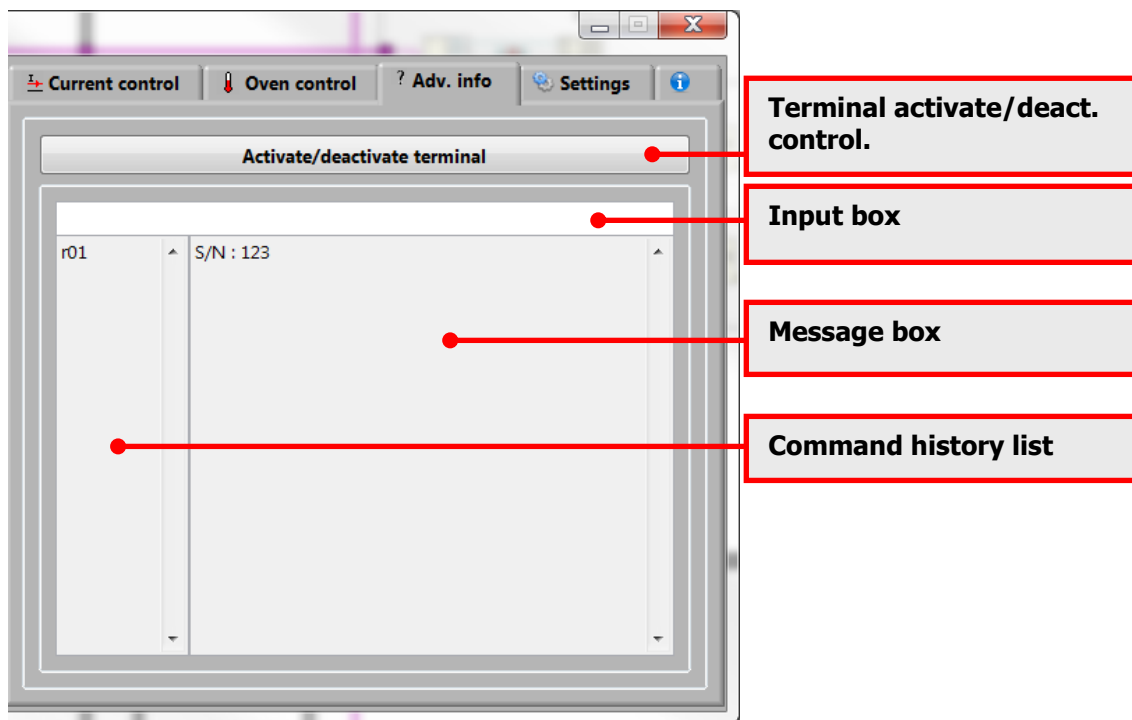


Figure 20: Advanced Info tab – terminal panel activated

The *terminal panel* consists of *input box*, *command history list* and *message box*.

#### Input Box

A command should be input in this box. Press Enter key when the input focus is inside this box to send out a non-empty command (empty command means spaces only).



In a terminal emulator, e.g. HyperTerminal, keystroke is sent out immediately. As a result, improper key sequence cannot be corrected. Unlike it, the command can always be revised inside the input box before sent out.

### Command History List

This box holds a list of recently used commands. Once a command is send through the input box, it will be appended to the end of the history list. To resend one of command, just double click with the left button of the mouse on the item.

### Message Box

Messages received from laser are displayed here. The content can be copied by using the windows hot key "Ctrl + C".

### 3.3.2.4 Settings Tab

The Settings tab consists of laser connection panel, hot keys panel and an info button (Figure 21), which shows the info dialog. Please refer section 4 for more information.

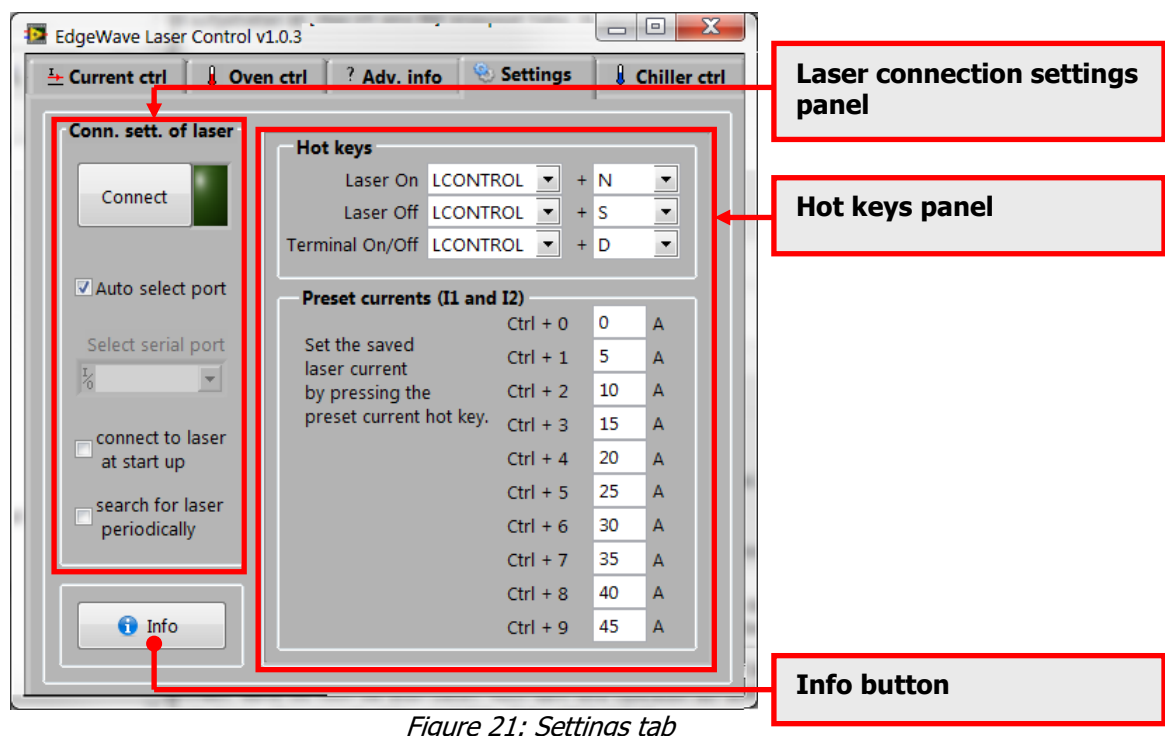


Figure 21: Settings tab

#### Laser Connection Settings Panel

Options for laser connection. Please refer section 4 for detailed information.

#### Hot Keys Panel

Configuration panel for hot keys. Please refer section 4 for detailed information.

### 3.3.2.5 Chiller Control Tab

The EdgeWave laser systems will be delivered in combination with a chiller. The most chiller have a serial port (RS232) to control it by a computer. In this tab (Figure 22) it is possible to control the chiller.

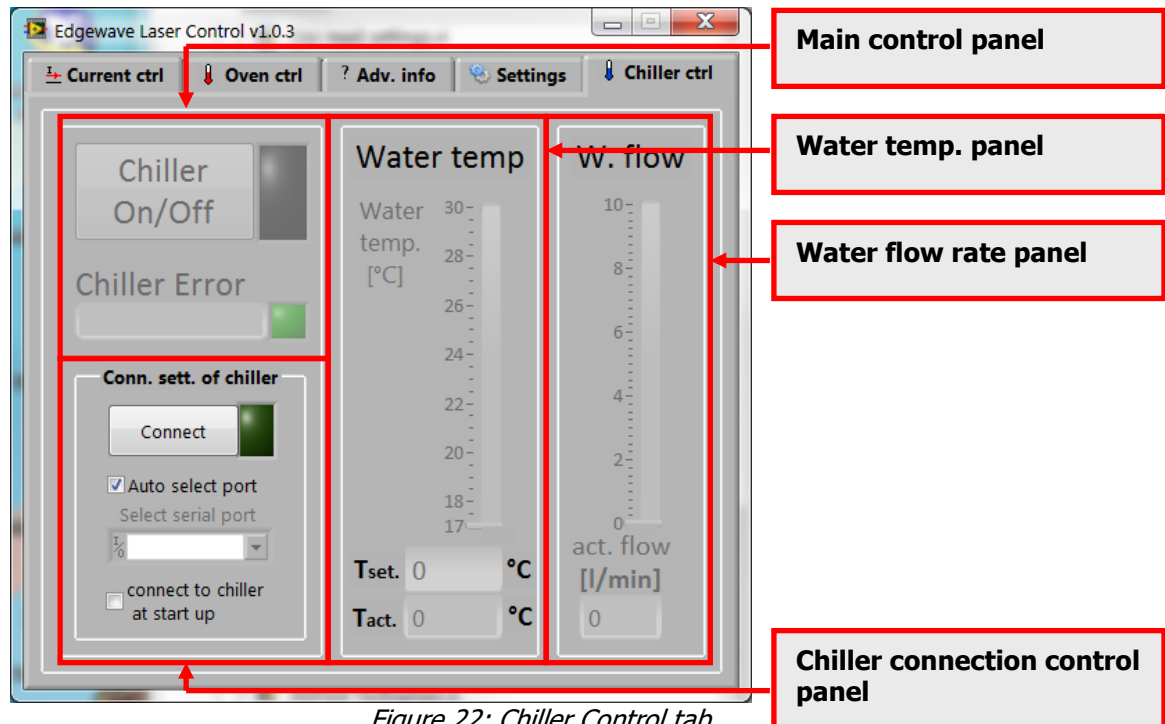


Figure 22: Chiller Control tab

#### Chiller Connection Control Panel

In this panel the connection settings can be changed and a manually connection to a chiller can be done. The settings will be saved at program end.

#### Connect Button

By clicking on *Connect* button the software will try to establish a connection to the chiller which is connected over a free serial port to computer.



The connection can be established only if the requirements in section 3.1 are fulfilled.

#### Auto Select Port Control

If enabled (checked) it detects the serial port automatically if a chiller is connected over a free serial port.

### Select Serial Port Control

Serial port number has to be specified correctly. If the *Auto select port* option is not checked. The software can enumerate the serial ports present in the computer which will be shown when click the down arrow in the drop-down list control.

### Connect To Chiller At Start Up

If enabled (checked) the software will try to connect to a chiller at program start up.

### Main Control Panel

It consists of a control element, a few indicators and message boxes.

### Chiller On/Off Control

The connected chiller can turn on or off by activating or deactivating this control.



If trying to turn off chiller and connected laser is on, a dialog will pop up to warn that the laser is on. It is possible to turn off laser and chiller at same time or cancel the operation.

### Chiller On/Off Indicator

It shows if chiller is on or off. The color of the indicator is green if chiller is on else black.

### Chiller Error box

On error or warning a message will be displayed here.



If the chiller is turned off, a message will be displayed and the *Chiller error* indicator will change its color to red.

### Chiller Error Indicator

On error or warning the color of the indicator is red, else green.

#### Water Temperature Panel

In this panel the cooling water temperature can be changed. *Tact.* and a slider show the actual outlet water temperature.

##### Tset

The temperature set point will read from chiller while connection and can be changed here. Nominal value is sent to the controller only after press Enter key in the edit box.



The water temperature can be set between 17°C and 30°C

##### Tact.

Tact. and a slider show the actual outlet water temperature

#### Water Flow Rate Panel

*Act. flow* and a slider shows the actual water flow rate in this panel.

## 4 Options

The Settings tab consists of laser connection, hot keys, preset panel and an info button (Figure 23), which shows the info dialog.

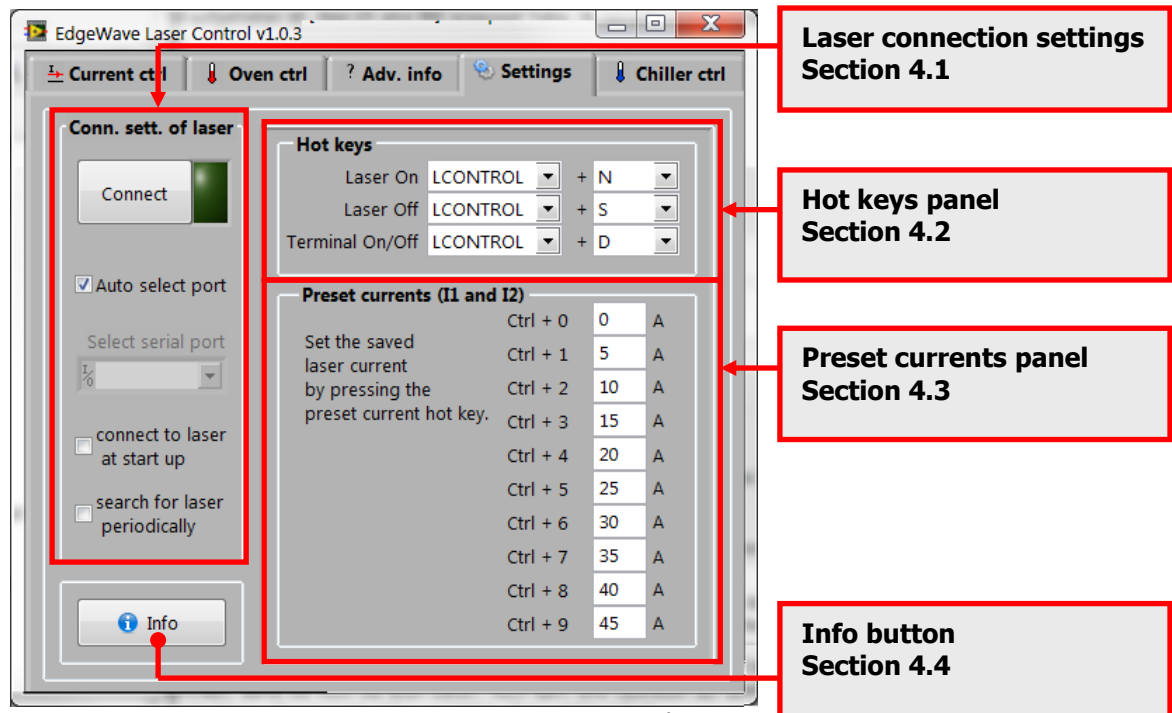


Figure 23: Settings tab

### 4.1 Laser Connection Settings

To establish a connection to laser, the following requirement must be fulfilled:

- An EdgeWave laser is connected to a working serial port of the computer.
- The RS232 cable used to connect the laser and the computer must be a 1-to-1 cable.
- Serial port number has to be specified in the laser control software correctly. If the *Auto select port* option is not checked.
- Emergency stop button of the laser has to be released.
- Safety loops on the rear panel of laser power supply is connected properly and the corresponding switches are closed. Please find more details about safety loops in the user manual of EdgeWave Laser.
- The Laser power supply is connected to the mains and switched on.



### Connect Button

To connect or disconnect a laser.

### Connection Indicator

If no laser is connected, the indicator is dark green. When the software tries to connect a laser, it changes its color to yellow. The indicator has the color bright green, if the software is connected to a laser.

### Auto Select Port

The Laser Control software enumerates the serial (RS232) ports present in the computer. If this option is activated, the software tries to open the ports and sends a command (r01 for serial number).

### Select Serial Port

The software has to know which serial port (RS232) is used for the connection to a laser.

The software can enumerate the serial ports present in the computer which will be shown when click the down arrow on the *select serial port* control.



The control element is deactivated (grayed out), if the *Auto select port* option is checked.

### Connect To Laser At Startup

When this option is enabled (checked), the software tries to connect laser at startup.

### Search For Laser Periodically

When this option is enabled (checked), the software tries to connect laser every five seconds until a laser connection can be established.



The software will not try to reestablish the connection if it is disconnected manually by the *Connect* button.

## 4.2 Hot Keys

Hotkey provides easier control in the software. To change the key combination used for certain function, use mouse scroll to select the key listed in the list box.

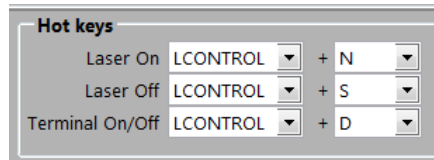


Figure 24: Hot keys panel



The software will NOT perform a conflicting check when assigning hotkeys. Please check all these fields carefully and avoid using the following hotkey which have been assigned to other functions:

Alt+X, Ctrl+O, Ctrl+D, Ctrl+C, Ctrl+V, F1 and Ctrl+0 ~ Ctrl+9.

### Laser On

Turn laser state to on by using the assigned hot key.

### Laser Off

Turn laser state to off.

### Terminal On/Off

Turn Terminal on or off.

## 4.3 Preset Currents

This feature provides a faster way to change the laser current easily. The most frequently used current values can be set here.

When Laser is connected, pressing Ctrl+0 ~ Ctrl+9 will set the nominal laser current to the value listed here. If the Laser has 2 branches, nominal currents of both branches (Current 1 and current 2) will be set according to the preset value simultaneously.

Preset currents (I1 and I2)		
Set the saved laser current by pressing the preset current hot key.	Ctrl + 0	0 A
	Ctrl + 1	5 A
	Ctrl + 2	10 A
	Ctrl + 3	15 A
	Ctrl + 4	20 A
	Ctrl + 5	25 A
	Ctrl + 6	30 A
	Ctrl + 7	35 A
	Ctrl + 8	40 A
	Ctrl + 9	45 A

Figure 25: Preset currents panel

### 4.4 Save Settings

The configuration of EdgeWave Laser Control is stored in the file called *Settings.ini* located at the same directory as the control software itself. The settings including the options introduced above will be loaded when launching Laser Control software. If *Settings.ini* cannot be found or corrupted, default settings will be loaded.

When the software is closed, all settings will be written into the setting file automatically. The software will create one if *Settings.ini* cannot be found and current Windows user has the privilege to create a file at the same folder.

Backup of the settings can be done by making a copy of *Settings.ini*.